

Peroxide Block

Description: Peroxide Block has been developed to use with immunolabeling techniques for the reduction of nonspecific background staining due to endogenous peroxidase.

Appearance: Clear liquid

Uses: Immunohistochemical procedures.

Limitations: Do not use past expiration date.

Availability:	<u>REF</u>	<u>Volume</u>
	ACA125	125 ml
	ACA500	500 ml
	ACA999	1000 ml

Storage: Store Peroxide Block at 2-8°C. Product is stable for 18 months from date of manufacture.

Procedure:

1. Incubate tissue section for 10 minutes at either room temperature or 37° C prior to application of the primary antibody. After incubation, rinse once in buffer.
2. *** For bulk staining, pour Peroxide Block in a covered staining tray and dip slides for 5 minutes. Replace with fresh reagent after 5-10 uses.

Precautions: Do not pipette reagent by mouth.
Avoid contact with skin and eyes.
Wash after use.
Observe all federal, state and local environmental regulations regarding disposal.

Storage: 2° C



8° C



ScyTek Laboratories, Inc.
205 South 600 West
Logan, UT 84321
435-755-9848
U.S.A.



Authorized Representative in Europe
(Regulatory affairs only)
EmergoEurope (31)(0) 70 345-8570
Molsstraat 15
2513 BH Hague, The Netherlands