

DAB AWAY

Contents:

Volume

125 ml DAB Away 1

125 ml DAB Away 2

250 ml Decolorizer

Storage: Store in dark at room temperature.

Procedure:

1. To prepare working solution, mix DAB Away 1, DAB Away 2 and DI water in the following ratio. Mix well (Note: solution will be dark purple).

80% DI water
10% DAB Away 1
10% DAB Away 2
2. Apply to stained glassware, staining pads, etc.
3. Allow to stand 5-10 minutes with occasional swirling.
4. Rinse with water.
5. Apply Decolorizer to surfaces for 1-2 minutes.
6. Rinse cleaned labware with water and allow to dry.

Precautions:

- * **All three components are acidic.**
- * Avoid contact with skin or eyes.
- * Wear protective gloves and eye covers.
- * In case of contact, wash area with copious amounts of water.