

Aqueous-Mount

Description: Aqueous-Mount is formulated for coverslipping sections stained with alcohol soluble chromogens, such as Fast-Red or AEC. No heating is required prior to use. Coverslips may be removed by soaking in water.

Availability:

<u>Item #</u>	<u>Volume</u>
AMT030	30 ml
AMT060	60 ml

Storage: Store at room temperature. Do not use past the expiration date.

Procedure: Apply several drops of Aqueous-Mount and coverslip. To remove coverslips, soak overnight in water.

Precautions: Avoid contact with skin and eyes. Wash after use. Observe all federal, state and local environmental regulations regarding disposal.