

DAB (Post) Enhancing Solution

Description: 3,3'Diaminobenzidine (DAB) is a widely used chromogen for immunohistochemical staining and immunoblotting. When in the presence of peroxidase enzyme, DAB produces a brown precipitate that is insoluble in alcohol. This solution has been developed to enhance DAB staining in a post staining procedure. The DAB will turn black after post-staining.

Availability/Contents:

<u>Item #</u>	<u>Volume</u>
DES015	15ml Dropper
DES125	125ml
DES500	500ml
DES999	1000ml

Storage: Room Temperature.

Procedure:

1. After DAB development, rinse in water.
2. Immerse tissue in DAB (Post) Enhancing Solution for 1-2 minutes.
3. Rinse tissue, counterstain, and coverslip.