

## DAB Enhancer Solution

**Description:** 3,3'Diaminobenzidine (DAB) is a widely used chromogen for immunohistochemical staining and immunoblotting. When in the presence of peroxidase enzyme, DAB produces a brown precipitate that is insoluble in alcohol. This solution has been developed to enhance DAB staining. The DAB will stain darker red/brown.

**Availability/Contents:**

<u>Item #</u>	<u>Volume</u>
ACM015	15ml Dropper
ACM125	125ml
ACM500	500ml
ACM999	1000ml

**Storage:** Room Temperature.

**Procedure:**

1. Prior to DAB incubation, add 20  $\mu$ l of DAB Enhancer to 1 ml of DAB ready-to-use reagent.
2. Stain with DAB reagent as usual.
3. Rinse tissue, counterstain, and coverslip.