

Hematoxylin, Gill's #2

Description: Gill's #2, hematoxylin is a stain with histological applications.

Results: Nuclei are blue.

Availability:

<u>Item #</u>	<u>Volume</u>
HG2500	500 ml
HG2999	1 liter
HG2010	10 liter

Procedure:

1. Stain nuclei 5-15 minutes with hematoxylin.
2. Place in Bluing Reagent (tap water will work in 2-3 minutes) until stain is blue.
3. Rinse in DI water.
4. Coverslip.