

## Hematoxylin, Gill's #3

**Description:** Gill's #3, hematoxylin is a stain with histological/cytological applications.

**Results:** Nuclei are blue.

**Availability:**

<u>Item #</u>	<u>Volume</u>
HG3500	500 ml
HG3999	1 liter
HG3010	10 liter

**Procedure:**

1. Stain nuclei 5-15 minutes with hematoxylin.
2. Place in Bluing Reagent (tap water will work in 2-3 minutes) until stain is blue.
3. Rinse in DI water.
4. Coverslip.