

## Mayer's Hematoxylin (Lillie's Modification)

**Description:** Mayer's Hematoxylin (Lillie's Modification) is a stain with histological applications.

**Results:** Nuclei are blue.

**Availability:**

<u>Item #</u>	<u>Volume</u>
HMM500	500 ml
HMM999	1 Liter
HMM010	10 Liters

**Procedure:**

1. Re-hydrate tissue as usual.
2. Stain nuclei 5 minutes with hematoxylin.
3. Place in Bluing Reagent (BRT500) until stain is blue.
4. Rinse in DI water.
5. Dehydrate, clear, and mount in permanent mounting medium.