

Alkaline Phosphatase Enhancer (Stable Solution)

Intended Use: To use as a signal enhancer in conjunction with alkaline phosphatase based immunostaining systems.

Storage: Store at room temperature and do not use beyond the expiration date stated in the bottle.

Procedure:

1. Once tissue sections have been incubated with alkaline phosphatase, wash them with buffer thoroughly.
2. Wipe the glass to remove excess buffer.
3. Incubate for 1 minute at room temperature.
4. Drain excess alkaline phosphatase enhancer and add substrate/chromogen without washing.

Precautions: Avoid contact with clothes and exposed skin. If accidentally contacted, flush with tap water immediately. Do not take internally.