

Bluing Reagent For Hematoxylin

Description: Bluing Reagent is designed to provide complete bluing of hematoxylin without damaging cell structure.

Results: Nuclei are blue.

Availability:

<u>Item #</u>	<u>Volume</u>
BRT999	1000ml
BRT500	500 ml

Procedure:

1. Stain with hematoxylin according to protocol.
2. Place in Bluing Reagent until nuclei are blue (1-2 minutes).
3. Rinse in DI water.
4. Dehydrate, clear, and mount in permanent mounting medium.